

KAMIYA BIOMEDICAL COMPANY

Elephant Serum Amyloid A (SAA) ELISA

For the quantitative determination of serum amyloid A (SAA) in elephant serum

Cat. No. KT-1905

For Research Use Only.

PRODUCT INFORMATION

Elephant Serum Amyloid A (SAA) ELISA Cat. No. KT-1905

PRODUCT

The **K-ASSAY®** Elephant Serum Amyloid A (SAA) ELISA is an enzyme immunoassay for the quantitative determination of serum amyloid A (SAA) in elephant serum. For research use only.

INTRODUCTION

Serum amyloid A (SAA) is a positive acute phase protein of approximately 12 kDa. Serum levels increase due to inflammation and infection in many species. In elephants, serum levels of 0 - 47.5 µg/mL have been reported for healthy elephants. Levels up to 410 µg/mL were found in sick elephants.

PRINCIPLE

The assay uses two elephant SAA peptide-specific antibodies. One is used for solid phase immobilization (microtiter wells). The other, conjugated to HRP, is used for detection. Diluted serum samples and calibrators are incubated in microtiter wells together with HRP conjugate for one hour. If present, SAA molecules are sandwiched between the immobilization and detection antibodies. The wells are then washed to remove unbound HRP conjugate. TMB is added and incubated for 20 minutes. If SAA is present, a blue color develops. Color development is stopped by addition of Stop Solution, changing the color to yellow. Absorbance is measured at 450 nm. The concentration of SAA is proportional to absorbance and is derived from a calibration curve.

COMPONENTS

- Anti-SAA coated plate (12 x 8-well strips)
- HRP Conjugate, 11 mL
- SAA stock
- 20X Wash Buffer, 50 mL
- Diluent, 2 x 50 mL
- TMB Reagent, 11 mL
- Stop Solution, 11 mL

MATERIALS REQUIRED BUT NOT PROVIDED

- Pipettors and tips
- Distilled or de-ionized water
- Polypropylene tubes
- Vortex mixer
- Absorbent paper or paper towels
- Plate incubator/shaker
- Plate washer
- Plate reader capable of measuring absorbance at 450 nm
- Graphing software

GENERAL INSTRUCTIONS

1. All reagents should be allowed to reach room temperature before use.
2. Reliable and reproducible results will be obtained when the assay is carried out with a complete understanding of the instructions and with adherence to good laboratory practice.
3. The wash procedure is critical. Insufficient washing will result in poor precision and falsely elevated absorbance readings.
4. Laboratory temperature will influence absorbance readings. Our ELISA kits are calibrated using shaking incubators set at 150 rpm and 25°C. Performance of the assay at lower temperatures will result in lower absorbance values

WASH SOLUTION PREPARATION

The wash solution is provided as a 20X stock. Prior to use, dilute the contents of the bottle (50 mL) with 950 mL of distilled or de-ionized water. Unused wash buffer may be stored at 4 °C for one week.

DILUENT

The diluent is provided ready to use. Do not substitute other buffers.

CALIBRATOR

1. The SAA stock consists of an elephant SAA polypeptide lyophilized in a stabilizing matrix. Prepare the 250 ng/mL calibrator as described on the vial label.
2. Label seven polypropylene tubes as 125, 62.5, 31.25, 15.63, 7.81, 3.91 and 1.95 ng/mL. Dispense 0.25 mL of diluent into each.
3. Pipette 0.25 mL of the 250 ng/mL SAA calibrator into the tube labeled 125 ng/mL and mix. This provides the 125 ng/mL SAA calibrator.
4. Similarly prepare the remaining calibrators by two-fold serial dilution. Use the calibrators within 30 minutes. The reconstituted 250 ng/mL stock should be stored at -20 °C for optimum stability.

SAMPLES

This kit was designed for measurement of SAA in serum. We found levels ranging from 5 µg/mL in healthy animals to >500 µg/mL in serum from sick elephants. We suggest testing each sample at dilutions of 250-fold and 10,000-fold to obtain at least one value within range of the calibration curve.

1. A 250-fold dilution can be obtained by mixing 2.0 µL of serum with 498 µL of diluent.
2. A 10,000-fold dilution can be obtained by mixing 12.5 µL of the 250-fold diluted sample with 487.5 µL of diluent.

PROCEDURE

1. Secure the desired number of 8-well strips in the cassette. Unused strips should be stored in a sealed bag with desiccant at 4 °C.
2. Dispense 100 µL of calibrators and samples into the wells. We recommend that calibrators and samples be run in duplicate.
3. Add 100 µL of HRP-conjugate to each well.
4. Incubate on a plate shaker at 150 rpm and 25 °C for one hour.
5. Empty and wash the microtiter wells 5x with 1x wash solution using a plate washer (400 µL/well).
6. Strike the wells sharply onto absorbent paper or paper towels to remove residual droplets if present.
7. Dispense 100 µL of TMB into each well.
8. Incubate on an orbital micro-plate shaker at 150 rpm at 25 °C for 20 minutes.
9. After 20-minutes, stop the reaction by adding 100 µL of Stop solution to each well.
10. Gently mix. It is important to make sure that all the blue color changes to yellow.
11. Read absorbance at 450 nm with a plate reader within 5 minutes.


RESULTS

1. Using graphing software, construct a calibration curve by plotting absorbance values of the calibrators versus log₁₀ of the SAA concentration.
2. Fit the calibration curve to a four-parameter logistic equation and derive concentrations for the samples.
3. Multiply the derived concentration of samples by the dilution factor to determine the concentration in the original sample.
4. If the A₄₅₀ values of samples fall outside the calibration curve, samples should be diluted appropriately and re-tested. Ideally, A₄₅₀ values of samples should fall within the mid-range of the calibration curve.

TYPICAL CALIBRATION CURVE


A typical calibration curve with absorbance at 450 nm on the Y-axis against SAA concentrations on the X-axis is shown below. This curve is for illustration only.

SAA (ng/mL)	A450
250	2.625
125	2.195
62.5	1.371
31.25	0.951
15.63	0.568
7.81	0.471
3.91	0.383
1.95	0.349

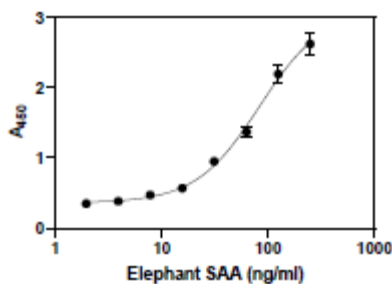

PERFORMANCE

Linearity: To assess the linearity of the assay, three samples containing SAA at concentrations of 11, 160 and 507 µg/mL were serially diluted with diluent to give values within range of the calibration curve.

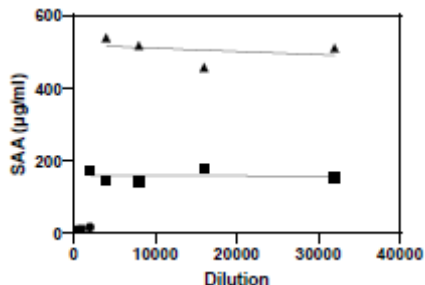

STORAGE

The kit should be stored at 4°C. The microtiter plate should be kept in a sealed bag with desiccant. Kits will remain stable until the expiration date.

FOR RESEARCH USE ONLY

KAMIYA BIOMEDICAL COMPANY

12790 Gateway Drive, Seattle, WA 98168
 Tel: (206) 575-8068 Fax: (206) 575-8094
 Email: LifeScience@k-assay.com
 www.k-assay.com