



KAMIYA BIOMEDICAL COMPANY

Biotin Labeling Kit - NH2

For the preparation of biotin-labeled antibody for enzyme immunoassay (EIA)

Cat. No. KT-221

For Research Use Only. Not for use in diagnostic procedures.

PRODUCT INFORMATION

Biotin Labeling Kit - NH2 Cat. No. KT-221

PRODUCT

Biotin Labeling Kit - NH2 is primarily used for the preparation of biotin-labeled antibody for enzyme immunoassay (EIA). NH2-Reactive Biotin, a component of this kit, has a succinimidyl ester group, and can easily make a covalent bond with an amino group of the target protein or other macromolecules without any activation process. Filtration Tube included in this kit is used for sample protein in removing small molecules such as Tris buffer and amine compounds that interfere with the assay or labeling reaction. The labeling process is very simple. Add the NH2-Reactive Biotin to protein solution on a filter membrane, and incubate at 37°C for 10 min. Excess biotin molecules can be removed by using a Filtration Tube. This kit contains all of the necessary reagents for labeling, including the storage buffer for conjugates.

COMPONENTS

- NH2-Reactive Biotin 3 tubes
- Reaction Buffer 500 µL
- WS Buffer 4 mL
- Filtration Tubes 3 tubes

Materials or equipment required but not provided

- Microtubes
- 10 µL, 200 µL adjustable pipettes
- Microcentrifuge
- DMSO
- Incubator (37°C)

CAPACITY

Three samples labeling

Sample requirement: Molecular weight >50,000; amount: 50-200 µg

PROCEDURE

1. Add 100 µL of WS Buffer and the sample solution containing 50-200 µg protein to a Filtration Tube. The volume of sample solution should be less than 100 µL. If the protein concentration is lower than 0.5 mg/mL, repeat Steps 1 and 2 until the total protein accumulation becomes 50-200 µg.
2. Pipette to mix and centrifuge at 8,000 x g for 10 min.
If the solution still remains on the membrane after the centrifugation, spin for another 5 min.
3. Add 10 µL DMSO to the NH2-Reactive Biotin, and dissolve with pipetting.
NH2-Reactive Biotin is on the bottom of the tube. Add 10 µL DMSO to the bottom of the tube, and pipette several times to dissolve.
4. Add 100 µL Reaction Buffer to the Filtration Tube, and then add 8 µL of NH2-Reactive Biotin solution to the Filtration Tube and pipette to mix.
If the amount of protein is 200 µg, add entire NH2-Reactive Biotin solution.
5. Incubate the tube at 37°C for 10 min.
6. Add 100 µL WS Buffer to the Filtration Tube, and centrifuge at 8,000 x g for 10 min. Discard the filtrate.
If the solution still remains on the membrane after the centrifugation, spin for another 5 min.
7. Add 200 µL WS Buffer to the Filtration Tube, and centrifuge at 8,000 x g for 10 min. Repeat this step one more time.
If the solution still remains on the membrane after the centrifugation, spin for another 5 min.
8. Add 200 µL WS Buffer, and pipette about 10 times to recover the conjugate. Transfer the solution to a microtube (not included in this kit), and store at 4°C.

We recommend using WS Buffer to recover the conjugate. You can choose any kinds of buffers appropriate for your experiment.

PRECAUTION

If the target protein solution contains other proteins with molecular weight larger than 10,000, such as serum albumin or gelatin, purify the protein solution, and use the purified target proteins for biotin labeling, because it might interfere the filtering or labeling reaction.

If the protein solution contains small insoluble materials, centrifuge the solution, and use the supernatant for the labeling.

The droplets which induced from the dry inhibitor of membrane, are occasionally found inside Filtration Tube while storing the kit at 4°C or after returning to room temperature. This phenomenon does not affect the performance.

STORAGE

Store at 4°C. This kit is stable for 1 year at 4°C before opening.

CAUTION: After a NH₂-Reactive Biotin is taken out from the seal bag, keep the unused NH₂-Reactive Biotin(s) in the bag, seal tightly and store at -20°C. Store the other components at 4°C.

Q & A

Q. Which protein can be labeled with biotin using this kit?

A. It can be possible to label protein that the molecular weight must be over 50,000 with reactive amino-groups.

Q. Can I use this kit to label antibodies which is commercially available?

A. Yes. However, if antibody solution contains other proteins such as serum albumin or gelatin, labeling reaction might be interfered by that protein. Purification of the antibody solution with affinity chromatography is necessary prior to use this kit. Contact us for the purification procedure, if you need.

Q. How long is the biotin-labeled protein stable?

A. Stability of conjugate depends on the protein itself. In the case of labeling for rabbit IgG, the labeled IgG is stable at 4°C for 2 months. However, for longer storage, add equal volume of glycerol to the sample solution and store at -20°C.

Q. How many biotin molecules are introduced to protein?

A. The number of conjugated biotin depends on the protein. In the case of rabbit IgG, 7 to 10 biotin molecules conjugate to each protein molecule.

Q. What is the minimum amount of protein that can be labeled using this kit?

A. We recommend using 50 µg as a minimum amount. Though 10 µg protein can be labeled using this kit, the background will be increased.

Q. Can I use this kit to label oligonucleotides or oligopeptides?

A. No. Oligonucleotides and oligopeptides may be too small to retain on the membrane filter of the filtration tube.

Q. Is there any notice for treatment of living cells with the Biotin-labeled protein?

A. We recommend using PBS including 2-10% FBS for preparation of cell suspension to maintain the best cell conditions.

Q. Does recovery buffer (WS Buffer) have harmful effect to living cells?

A. No. WS Buffer contains stabilizing agent (surfactant) that is controlled of its concentration without cytotoxicity. If you are concerned about the additive in WS Buffer, you can use your own buffer currently used instead of WS Buffer.

FOR RESEARCH USE ONLY

KAMIYA BIOMEDICAL COMPANY

12779 Gateway Drive, Seattle WA 98168

Tel: (206) 575-8068 Fax: (206) 575-8094

Email: LifeScience@k-assay.com

www.k-assay.com